

Laptop Mount Assembly Instructions

Page 1


The Laptop Mount allows you to use a Gazepoint eye tracker with a laptop.

Laptop Mount Kit includes:

Laptop Mount


Black M4 Thumbscrews (2)

We recommend that you first install the Gazepoint installer software before mounting the eye tracker.

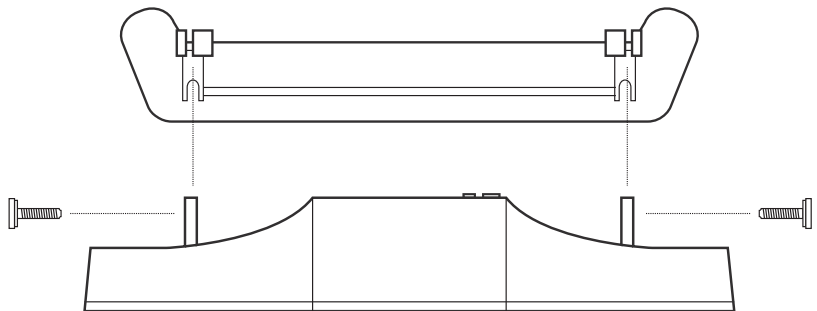

1

Position the Laptop Mount so that the flared ends are to the back of the eye tracker. Push the eye tracker mount tabs into the slots of the Laptop Mount and thread the M4 Thumbscrews to tighten.

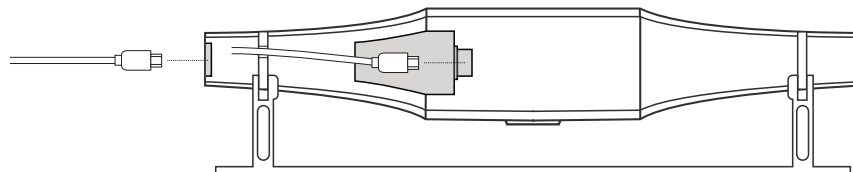


Gazepoint Research Inc.
418 – 2912 West Broadway
Vancouver B.C.
V6K 0E9 Canada

T +1 408 883 4293 (GAZE)
E info@gazept.com
W www.gazept.com

2

Connect the USB and power cables to the eye tracker and place on the keyboard as close to the lower edge of the screen as possible and angle upwards to approximately 45° towards your face and tighten M4 Thumbscrews.

- 3 Plug in the USB power and data cables to your computer.

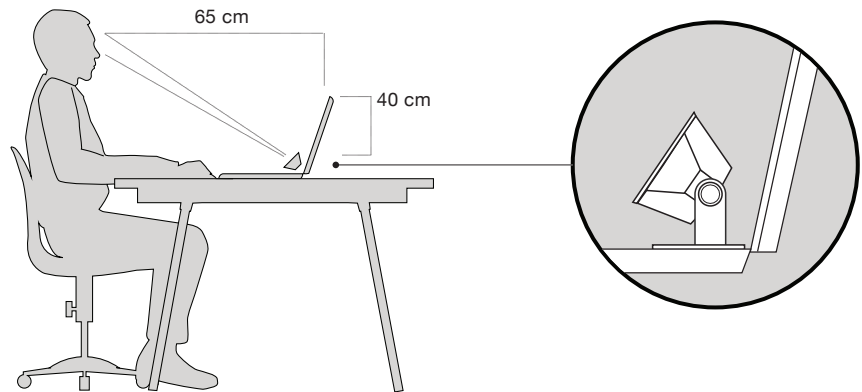

- 4 Position Laptop so that the eye tracker is approximately 65cm (25") from user.

Gazept Research Inc.
418 – 2912 West Broadway
Vancouver B.C.
V6K 0E9 Canada

T +1 408 883 4293 (GAZE)
E info@gazept.com
W www.gazept.com